

R.T.U X-Gal IPTG Solution (1000X)

Cat No: PT-BL006

Size: 2mL

Concentration: 125 mg/ml X-Gal, 119mg/ml IPTG

Description:

Protech PT-BL006 X-Gal IPTG Solution (1000X) is a ready-to-use solution for blue/white selection of cloning constructs in standard lacZ systems. X-Gal IPTG Solution (1000X) is composed of 125 mg/ml X-Gal and 119 mg/ml IPTG in a non-hazardous and non-toxic solvent.

General Protocol:

1. Cool agar to about 50°C.
2. Add 100 µL R.T.U. X-gal IPTG Solution (1000X) to every 100 mL of agar. Mix it gently.
3. Pour agar into sterile petri dishes.
4. Allow agar to solidify.
5. Plates could be used immediately or stored at 2 - 8°C and protected from light.

Storage:

Stable for a year when stored at -20°C.

*****Research Use Only*****

Please do not hesitate to contact us if you have any questions.

Manufactured for and distributed by Protech Technology Enterprise Co., Ltd.

TEL : +886-2-2655-7677 / FAX : +886-2-2655-7601 / Toll Free : 0800-231-530

E-MAIL: service@bio-protech.com.tw ; tech@bio-protech.com.tw

WEB: <http://www.bio-protech.com.tw>