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Free!

MMLV Reverse Transcription kit 100 reactions 特價 \$4,800 (20%off)

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
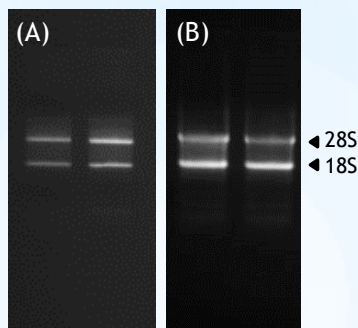
High-quality total RNA extracted from (A) 1.2×10^6 human epidermis cells and (B) rat tissue

另有miRNA純化套組: **SanPrep Column microRNA Miniprep Kit** (BioBasic Cat. No. SK8811) for the isolation of miRNA, siRNA and snRNA (<200 nt).

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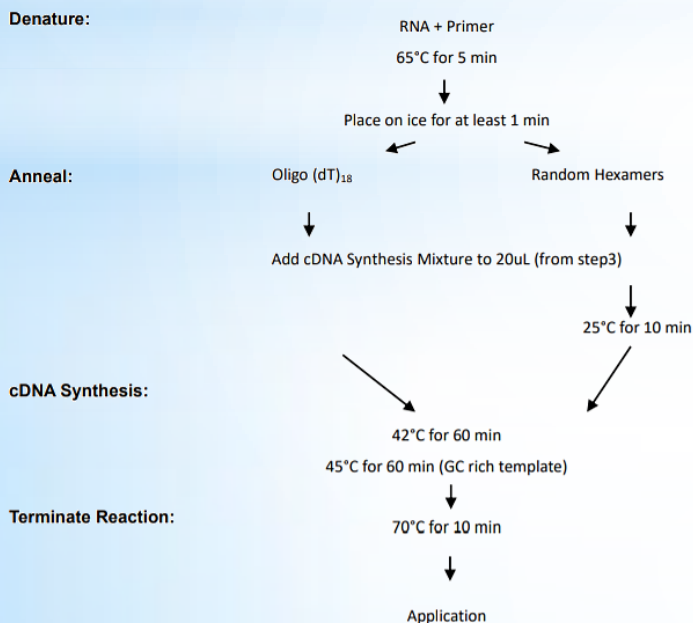MMLV Reverse Transcription Kit Flow Chart

Gene-Spin™ Total RNA Kit Flow Chart

	Animal	Plant	Cell	Bacteria
Homogenize Sample	20-30 mg	100 mg	5×10^6 $\sim 1 \times 10^7$	$< 1 \times 10^9$ 100 μL Lysozyme RT, 3-10 min
RNA Lysis/2ME Solution	600 μL	450 μL 60 °C, 3 min	600 μL	350 μL
Spin Lysates	14000 x g 3 min	14000 x g 3 min	-	14000 x g 2 min
Lysates + EtOH Mix Well	70% EtOH 600 μL	100% EtOH 225 μL	70% EtOH 600 μL	100% EtOH 250 μL
Bind RNA	Load sample 14000 x g 1 min			
Wash I	500 μL 14000 x g 1 min			
DNase I Digestion	2 μL DNaseI (2U/μL) 80 μL DNaseI Buffer RT, 15 min			
Wash I	500 μL 14000 x g 1 min			
Wash II	1st. 600 μL 14000 x g 1 min			
	2nd. 600 μL 14000 x g 1 min			
Remove Residual EtOH	14000 x g 3 min			
Elute RNase-free H ₂ O	30~50 μL 14000 x g 1 min			

MMLV Reverse Transcription Kit (PT-RT-KIT)

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
Order Information


Product Name	Size	Contents	Cat. Number
Gene-Spin Total RNA Purification Kit	50 reactions	Spin column, RNA lysis solution, DNase I enzyme and buffer, RNA wash solution I & II, nuclease-free water	PT-RNA-MS-50
MMLV Reverse Transcription kit	100 reactions	MMLV RT enzyme, buffer, dNTP, oligo dT ₁₈ primer, random hexamers, DEPC water and RNase Inhibitor	PT-RT-KIT

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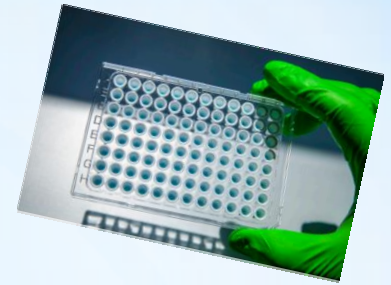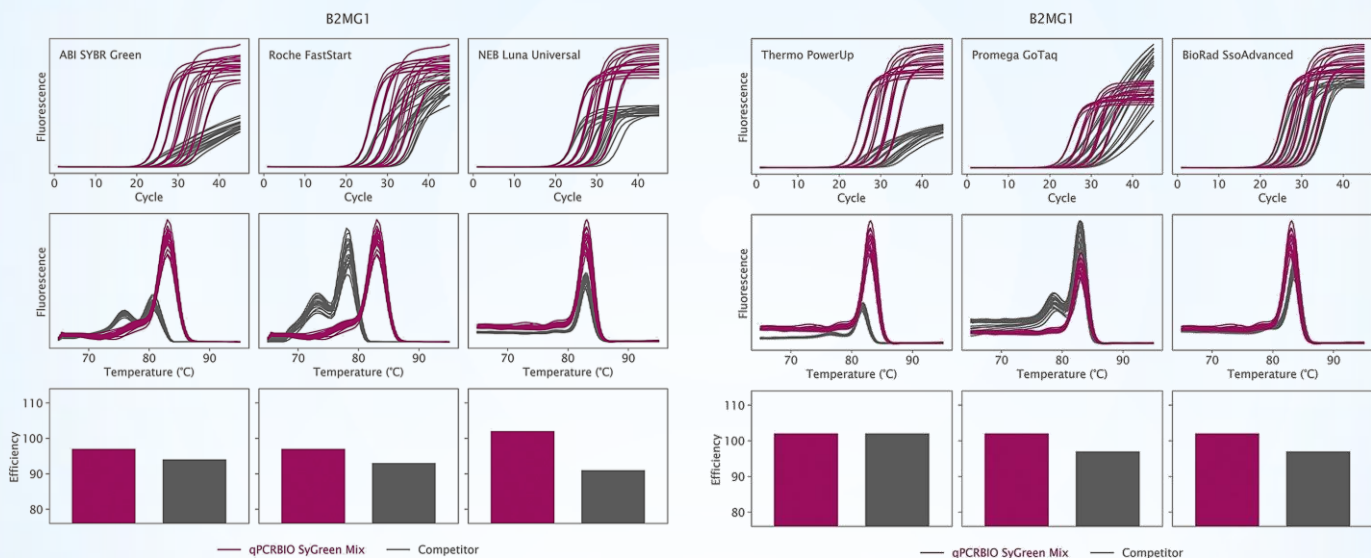Comparison of qPCR BIO SyGreen Mix (purple curves) with six other competitors. Using mouse cDNA as template, beta-2 microglobulin was amplified according to conditions set by each of the competitors. From top to bottom panels were: amplification curves, melting curves and amplification efficiencies.

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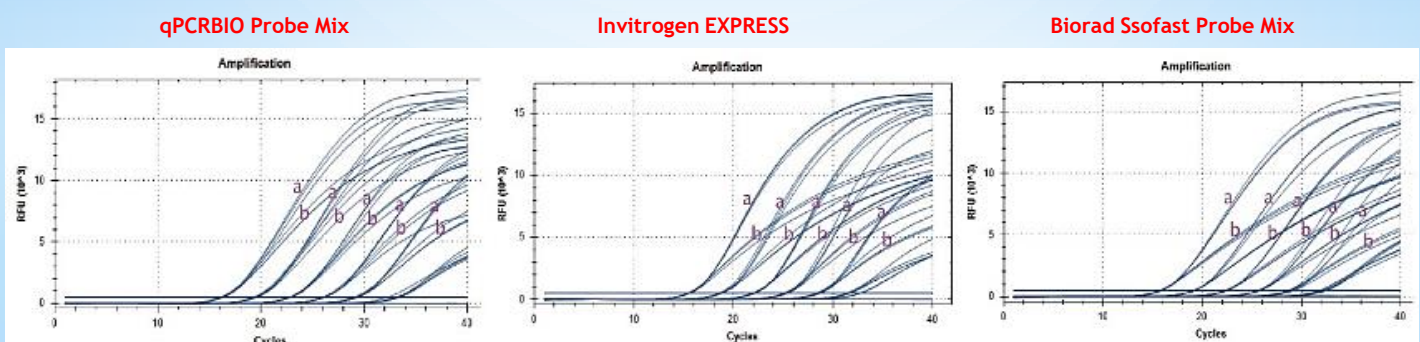
qPCR BIO Probe Mix has the least multiplex inhibition effect!

TaqMan probe amplification of human LIMK1 gene in a=singlex or b=quadplex (ACVR2B, LIMK1, ACVR1B and CDK7). Amplification is less inhibited in multiplex with qPCR BIO Probe Mix comparing with the other two mixes.

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