

Product No. 25969

Phenol, Saturated with TE Buffer

Components

Volume	Components	
	Phenol, Saturated with TE Buffer	Tris buffer
100 ml (Product No.25969-54)	100 ml	6.5 ml
400 ml (Product No.25969-96)	400 ml	26 ml

Protocol

High pH(pH 7.9±0.2)

1. Add Tris buffer into the product bottle. Shake to mix gently.
2. Leave it for 2-4 hours until two phase separations. [pH of lower phase (Saturated phenol) is adjusted to 7.9±0.2]
3. Use the lower phase. Be careful not to mix upper phase.

Low pH(pH 6.6±0.2)

1. Not add Tris buffer.
2. Use lower phase after leaving until two phase separations. Be careful not to mix upper phase.
[pH of lower phase (Saturated phenol) is adjusted to 6.6±0.2]

Storage

Refrigerator