

Date: 2016.08
Version: 05_04

4X Protein Loading Dye

Catalog number : SA-PRO-4X-DYE	Size : 100 ul
PT-PRO-4X-DYE	1 ml
PT-PRO-4X-DYE-50ML	50 ml

Components : (For 4X concentration)

200mM Tris-HCl (pH6.8), 4% 0.57M 2-mercaptoethanol, 8% SDS, 0.04% Bromophenol blue, 40% Glycerol

Description :

Incubate Protein sample at > 95°C for 2-5 minutes in 1X loading dye before it electrophoreses through SDS-PAGE gel.

Storage condition :

The solutions are stable at -20°C. However, for long term storage, we recommend storing in aliquots at -70°C.

For Research Using Only.

Please do not hesitate to contact us if you have any questions.

Manufactured for and distributed by Protech Technology Enterprise Co.,Ltd

TEL : +886-2-2655-7677 / FAX : +886-2-2655-7601 / Toll Free :
0800-231-530

E-MAIL: service@bio-protech.com.tw ; tech@bio-protech.com.tw

WEB: <http://www.bio-protech.com.tw>