波仕特Western Blot專刊

1

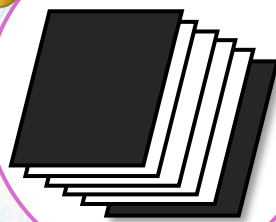
Sample Preparation

2


Gel Electrophoresis

3

Membrane transfer

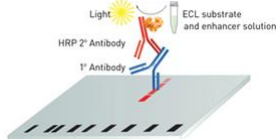4

Immunodetection

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
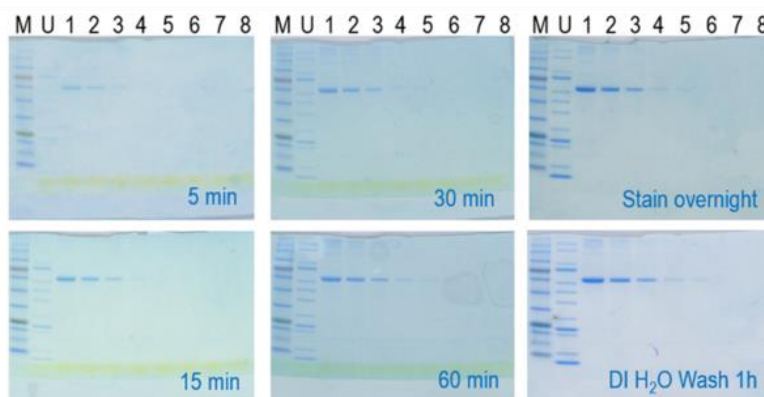
M: Prestain ladder, U: Unstain ladder, Lane1: BSA 2000 ng, Lane2: BSA 1000 ng, Lane3: BSA 500 ng, Lane4: BSA 100 ng, Lane5: BSA 50 ng, Lane6: BSA 10 ng, Lane7: BSA 5 ng, Lane8: BSA 1 ng

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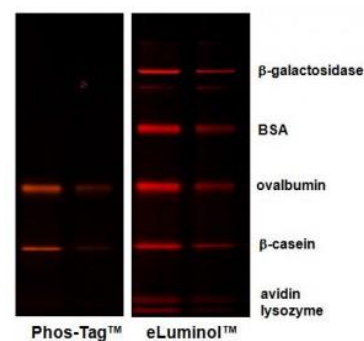Condition:

Electrophoresis gel Bullet PAGE One Precast Gel 5-15% (#13079)
 Samples
 Protein Ladder One (#09547), 5 µl added to each lane
 Protein Markers (10x) (#29458), 2 µl added to each lane
 Protein Markers (10x) (#29458), 3x diluted, 2 µl added to each lane
 Protein Markers (10x) (#29458), 10x diluted, 2 µl added to each lane
 Protein Markers (10x) (#29458), 30x diluted, 2 µl added to each lane
 Membrane PVDF Transfer conditions: 10V (constant voltage), 10 min; Detection CBB Stain One (#04543)

Transfer buffer	Bullet Semi-Dry Transfer One	Kyhse-Andersen	Towbin
Conditions	10V (constant), 10 min		
Membrane image			

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Sample: Hela cell extract; (left) 2 µg, (right) 0.4 µg
 Blocking: Incubated for 60 minutes with 5% skim milk
 Primary antibody: β-Actin antibody manufactured by Santa Cruz, #sc-47778, X 1,000 dilution
 Secondary antibody: Goat anti-mouse IgG(H+L) secondary antibody manufactured by Novous, NB7574, X 40,000 dilution
 Detection: Chemi-Lumi One Super manufactured by Nacalai, #02230, 5-minute exposure time

Signal Enhancer HIKARI for Western Blotting and ELISA

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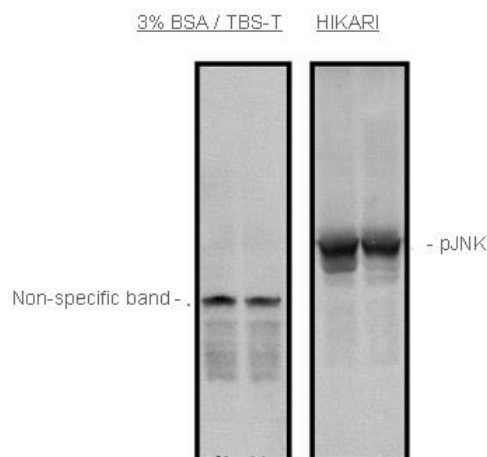
Primary Ab: Anti-Rac1 (1: 1,000)
Secondary Ab: Anti-Mouse IgG-HRP (1: 2,500)
Sample: Rat cortical primary neuron
Detection: ECL

Data courtesy of Dr. K. Ishizuka,
Department of Psychiatry, Johns Hopkins University School of Medicine

Primary Ab: Anti-pJNK (1: 1,000)
Secondary Ab: Anti-rabbit IgG - HRP (1: 5,000)
Sample: Mouse Embryonic Fibroblast (MEF)
Detection: SuperSignal West Pico

Data Courtesy of Dr. S. Matsuzawa,
Signal Transduction, NCI Cancer Center, Burnham Institute for Medical Research

Bovine Serum Albumin

Appearance
White to off-white coloured, lyophilized powder.

- Ultra High Purity 超高純度
- Low Immunoglobulin 低免疫球蛋白
- Low Endotoxin 低內毒素
- Low Fatty acid content 低脂肪酸含量
- Protease free classified 無蛋白酶
- Bovine virus free 無牛病毒
- Consistent batch to batch properties 批次品質穩定
- 詳盡的COA
- 完整的產地證明COO

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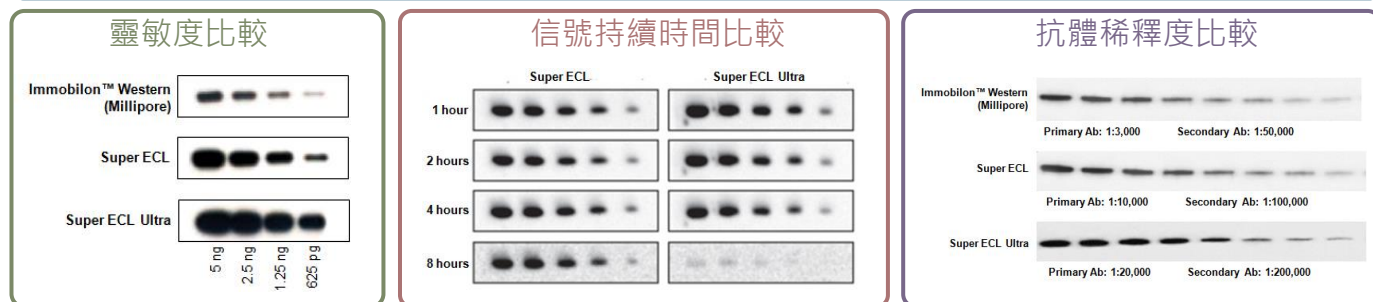
Experiment conditions:

Sample: Hela cell extraction, 2 times dilution started from 3 ug; 1 Ab: Anti-GAPDH manufactured by Novous, #NB300-322, X 5,000 dilution; 2Ab : Anti-rabbit IgG, HRP-linked whole Ab Donkey manufactured by Cytiva, #NA934, X 10,000 dilution. Detector for chemiluminescence : LAS-3000 Super mode

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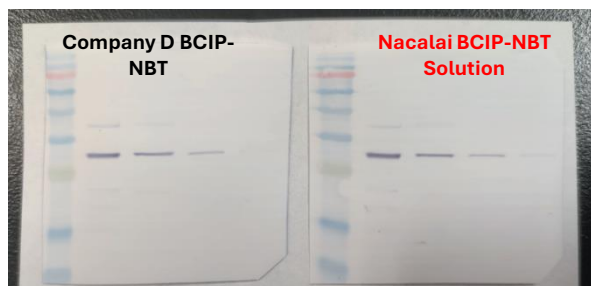
Western blotting 檢測

Experiment conditions

Sample : Protein from mouse brown adipose tissue (From left 30 µg, 10 µg, 3 µg, 1 µg)
Blotting : 5% skim milk/tTBS at room temperature for 1 hour.
Primary antibody : Anti-UCP1 antibody (Rabbit) (Self-preparation) 2,000x diluted
Diluted using 5% skim milk/tTBS, refrigerated overnight.
Secondary antibody : Anti-Rabbit IgG (Fc), AP Conjugate(Promega #S373B) 5,000x diluted
Diluted using 5% skim milk/tTBS, left at room temperature for 1 hour.
Detection : 1 minute for each staining solution.

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產品名	原廠編號	SIZE	說明
Super ECL Western Blotting Detection Kit	FP300 & FP301	100/500mL	HRP冷光Substrate，靈敏度高，信號持續時間長
Super ECL Ultra Western Blotting Detection Kit	FP302 & FP303	100/200mL	HRP冷光Substrate，超靈敏
BCIP-NBT Solution	19880-84	100 mL	即用型AP 受質溶液，可用於WB和組織細胞呈色(藍紫色)
TMB Solution for Western blot	18186-24	200 mL	即用型HRP-based western blot呈色劑